

## **OOSAFE® SURFACE DISINFECTANT**

Safe disinfection and cleaning for laboratory surfaces and floor. It is also compatible with glass, metal, rubber, plastic and wood.



- Economical: I:100 water dilution is required while many other disinfectans require I:10 water dilution.
- MEA Tested: Only disinfectant which comes with MEA Test Result for each lot.
- No VOC: No alcohol, no VOC release.
- Long Shelf Life: 3 years after manufacturing.
- No Waste: Can be used till last drop until expiry date.
- Cleaning and Disinfection: As a result of non-ionic soap content, it can clean and disinfect at the same time. Suitable for use in operating rooms since it can dissolve dried blood stains.
- Fast Disinfection: Complete disinfection after I5 minutes of application.
- Scientists Approved: It has been used by leading IVF clinics in more than 65 countries.
- Quality: Each bottle has security seal which assures safe transportation and quality controlled product. 2 liter disinfectant is packed in individual box together with dose cup.
- Microbiological Efficacy: Bactericide, fungicide (Candida), algicide and selectively virucide.
- Usage: Can be used even in laboratories which do not have enough ventilation, no VOC release after application.
- Application: Dilute Oosafe Disinfectant I:100 with water before use and wait I5 minutes after application. If it is used for laboratory floor, please remember to rinse the floor I5 minutes after application, otherwise it may be slippery.
- Active Ingredients: Quaternary ammonium compound (propionate):
  6.3 mg/g



